Timer setting “RODAC-WEIGHT” – Code n. 15496 INSTRUCTIONS MANUAL

International PBI S.p.A. Via Novara, 89 – 20153 Milano – Italy – Tel + 39 02 48779-1 – Fax + 39 02 40090010
E-mail: info@internationalpbi.it – www.internationalpbi.it

INSTRUCTIONS MANUAL

READ CAREFULLY THIS HANDBOOK

SAFETY INSTRUCTIONS

- Use the device just for the indicated purposes.
- The Device must be correctly installed according to manual instructions, before starting any operation.
- Never disconnect the device pulling electrical cable from the socket; .
- Replace immediately electrical cable when damaged ; never use a damaged or worn electrical cable.
- On removing the item connected to the electrical system pay attention to the electrical connection cable
- Always disconnect from electrical system on:
 - A. Repairing or maintenance; these operations must be carried out by qualified staff;
 - B. Keep the unit clean;
 - C. Use original spare parts and accessories for any replacement;
- Do not use this device in the presence of gaseous explosive.

INSTALLATION GENERAL CAUTION

CAUTION

When the electrical system socket is not compatible with device socket, replace it paying attention to codified cable colours.

CAUTION

Be aware that the voltage and frequency of the electrical system are available with device electrical income.

The device must be connected to a ground wire. For any question please contact an electrician.

CAUTION

Before starting-up the device be sure that ventilation grids are totally free to operate and not covered by any barriers.

INDEX

1. Description	pg. 4
2. Time setting	pg. 4
3. Sliding of timer	pg. 4
4. Sliding of a contact plate	pg. 4
5. Sterilization	pg. 4

1. DESCRIPTION

MINUTES
LCD DIGITAL DISPLAY
SECONDS

MINUTES BUTTON
SECOND BUTTON
START/STOP BUTTON

Back
BATTERY COVER
MAGNET

2. TIME SETTING

10 seconds contact time setting

1. Press the “second” button and keep it depressed until reaching the figure 10 on the digital display.
2. Push the start/stop button once to begin the countdown.
3. When the timer counts down to “00” a beeping alarm will sound. It will continue for 30 seconds or until the timer is turned off by pressing start/stop button.
4. The setting of 10 seconds is memorized and the cycles can be repeated as often as desired.
5. The timer uses a 1,5 G-13 battery.

3. SLIDING OF TIMER

The timer must be pressed in its rail-housing and fixed (if necessary) by the fixing side screw.

4. SLIDING OF A CONTACT PLATE

The contact plate, complete of lid, must be slided on the other side of the “Rodac-Weight”.
The position of fixing pins can be adjusted for contact plate of different producers.

5. STERILIZATION

The “Rodac-Weight” may be sterilized by autoclaving (the timer is not autoclavable!).

**To the attention of: Technical department International pbi S.p.A.
Via Novara, 89 - 20153 - Milano - Tel + 39 02/48.779.1-Fax +39 02/400.900.10**

AUTHORISATION RETURN NUMBER TO BE ASKED TO THE EXPORT DPT

Company:		
Address:		
City:	Zip Code	State
Phone	Fax	Email
Name to contact:		
Lab/Division:		
<input type="checkbox"/> return	<input type="checkbox"/> request of technical service	
Equipment:		
Model	Serial Number:	
for <input type="checkbox"/> calibration	<input type="checkbox"/> repair	<input type="checkbox"/> other (specify)
NOTE OF MALFUNCTIONING		
<input type="checkbox"/> ELECTRICAL PART		
<input type="checkbox"/> ELECTRONICAL PART		
<input type="checkbox"/> MECHANICAL PART		
<input type="checkbox"/> HYDRAULICS PART		
NOTES:		
ATTENTION: THE INSTRUMENT MUST BE SENT TO INT.PBI C/O STATIM VIA SAN GIUSTO 85 20153 MILANO ITALY		
SIGNATURE	NAME	DATE

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WARRANTY

WARRANTY CONDITIONS

International PBI warrants this product to be free of defects in materials and functionality
The repairs will be effected by International PBI or by authorized dealers as follows:

- Faulty components free replacement
The labour is always at customer's expense
- The equipment will be sent at buyer's expenses and risk and the return will be ex-works
- This warranty does not apply to damages resulting from tamperings, abnormal use and installation and does not cover transport damages
- International PBI shall not be liable for any commercial damages, whether incidental, consequential or otherwise derived from the use of instruments

ATTENTION

TO OUR NEW CUSTOMER

As a user of our equipment you are covered for a 12 months guarantee for all spare parts as a result of mechanical failure of the machine during normal use. In order to provide you with first class service, it is essential for us to know all your references and which equipment you have. We kindly ask you therefore to fill in the attached guarantee and return the card to us by first class mail.

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International pbi S.p.A.
Via Novara, 89 - 20153 Milano

HEREBY DECLARES

under its sole responsibility that the following product

Timer setting “Rodac-Weight” Code: 15496

- ◆ the following harmonized standards have been applied
EN 61010-1/2001, EN 61326-1/2000

- ◆ EC Directive 2002/95/EC of the European Parliament: “Restriction of the use of certain hazardous substances” (RoHS) for electric and electronic devices

The EC RoHS Directive has been released for the restriction of the use of certain hazardous substances in electric and electronic devices and basically says that with July 1st 2006 new productions of electric and electronic devices may not contain lead, mercury, cadmium, hexavalent chromium, polybrominated biphenyl (PBB), respectively polybrominated diphenylether (PBDE).

As a consequence of this EC Directive we have to introduce unleaded technology in our production and we have to make adaptations concerning the use of lead, mercury, cadmium, hexavalent chromium, PBB, PBDE according to the defined permissible values. The adaptation of our products to the RoHS conformity is confirmed.

- ◆ EC Directive 2002/96/EC of the European Parliament: “Disposal of old products”

Our products are designed and manufactured with high quality materials and components which can be recycled and reused according to the EC Directive 2002/96/EC. The user should inform himself about the separate collection system for electrical and electronic products. He should act according to the local rules and do not dispose of the old products with the normal household waste. The correct disposal of old products will help to prevent potential negative consequences for the environment and human health.

- ◆ and conforms with the technical requirements of the following directives
Machines directive 98/37/CEE; Low voltage directive 73/23/CEE
Electromagnetic compatibility directive 89/336/CEE

Il Presidente
Roberto Ligugnana

Milano 2008